

How To Fill A Pipet

1. Wash the pipet with water
2. Rinse the pipet with the solution that's going in the pipet
3. Wipe the outside with a wipe
4. Using a bulb (you don't push the bulb on to the pipet. Just gently press them together), suck the liquid to just past the mark - Don't get solution in the bulb. This is very bad !
4. Gently rest the tip of the pipet on the bottom of the flask. Remove the bulb, Using a dry index finger (not, I repeat not, your thumb), let the liquid out until the meniscus lies on the mark
5. Wipe the outside of the pipet
6. Run out the liquid down the side of the flask
7. When the pipet is 'empty', count to three and the gently touch the tip of the pipet to the surface of the liquid.